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TPP cryo tubes are made of polypropylene (PP) with external thread and star feet are designed for storage of biological material and samples down to -196°C . The screw caps with external thread reduce the risk of contamination and cross-contamination.

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The cryo tube is intended for single use.

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Note:

Follow national regulations when handling biological material, use appropriate protective clothing

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Take note of the regulations of aseptic working during the work process.

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TPP recommends storing the tubes exclusively in the gas phase of nitrogen (LN_2). If stored in the liquid phase, LN_2 could penetrate. During thawing process, the rapid evaporation would then cause a build-up of internal pressure. This could result in an explosion of the tube and release of biologically hazardous substances.

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Instructions:

- *Check the packaging material and the product for intactness. Use only faultless material.*
- *Check the expiry date (EXP) on label and packaging. Only use products with valid EX.*
- *Open the packaging in a sterile environment and remove a complete system for use.*
- *For a firm and safe stand during the work, place the cryotubes in the appropriate rack (#99016).*
- *Remove the cap and fill according to your laboratory routine. Please refer to the recommended 80 % and maximal 90 % filling quantity to allow the sample volume to expand.*

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Storage in Minus Temperatures:

- *When storing the tube in the minus temperature range, first test the multitude of influencing factors under routine conditions.*
- *Temperatures below 0 °C result in reduced elasticity for polypropylene. Do not stress frozen tubes mechanically such as dropping or hitting. Periodically test under routine conditions for potential issues that could damage your samples.*
- *Only use "freezing containers" for controlled freezing of cells before converting the cryotubes into the gas phase of a nitrogen tank.*

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Advice and Recommendations:

- *The graduation serves as a reference for the filling quantity and is not an absolute value.*
- *A labeling area and colored cap inserts allow identification of samples.*
- *A detailed freezing protocol for cells can be found on our homepage www.tpp.ch*

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Accessories:

- *Cryo Box L, # 99014 / Cryo Box S, # 99015*
- *Cryo Rack # 99016: The star foot in combination with the TPP rack allows one-handed operation.*
- *Colored cap inserts for identification*

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Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch.

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Kryorack # 99016

Instructions:

Designed for storage of cryo tubes and most other similarly designed vials in ultralow temperature.

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